

Kimtech™ A5 Sterile Boot Cover with Grasp Ties

Donning Technique

1

Select your boot cover size; inspect package for vacuum seal; tear along notches toward the product to open; flex package to separate boot covers.



2

Pull out one boot cover by the cuff.



3

Grasp ties with one hand and make sure they don't touch the floor.



4

Open up boot cover; point toe toward opening and pull up over calf. Hold ties to prevent touching the floor.



5

Hold up foot or rest on bench. Grasp ties at front of boot cover and wrap over shoe/boot; then wrap ties behind ankle and tie in front.



6

Place donned boot on the clean side.



7

Pull the remaining boot cover out of package by the cuff and repeat steps 3, 4 and 5. Proceed with final gowning steps.

