

Kimtech™ A5 Sterile Cleanroom Apparel with Clean-Don™ Technology: Gowning Technique

KIMTECH™

BEFORE GOWNING

- 1 (Pre-Entry)** Don Hair Net, Shoe Covers (and Beard Cover if applicable) after removing all jewelry and cosmetics.
- 2 (Gowning)** Wash hands and don first pair of sterile gloves. Sanitize gloves after donning each article if required.
- 3** Don mask and hood assuring a secure fit. (Integrated hood and mask incorporates both components into one.)

BEGIN GOWNING

Snaps allow gathered-up arms and legs to expand during gowning.



After the Kimtech™ A5 Sterile Cleanroom Apparel is gowned, add boot covers.

Complete gowning by adding goggles and a second pair of sterile gloves.



Visit our website at www.kimtech.eu
Technical help: kimtech@kcc.com