


Kimtech™ A5 Sterile Integrated Hood and Mask

Donning Technique



1

Select your integrated hood and mask; inspect package for vacuum seal; tear along notches toward the product to open.

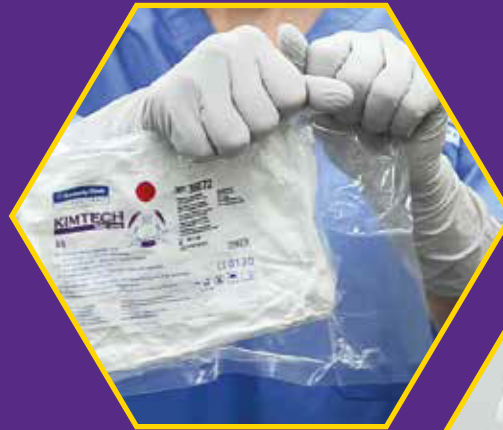

2

Pull out integrated hood and unfold to locate the blue signal indicator line.



3

Centre your chin just above the blue line then roll over your forehead.



4

Grab ties with both hands then extend arms out in a circular motion to position under arms.



5

Pull ties across chest and secure as needed.



6

Adjust front flaps to comfortable position, reach up underneath the bib of the hood, touching sterile, and pinch mask wire over the bridge of nose to ensure a secure fit.

